



Normal Mouse
serum blocking buffer

NB-23-00187-1

NB-23-00187-2

Normal Mouse serum blocking buffer

#Cat: NB-23-00187-1

Size: 100ml

#Cat: NB-23-00187-2

Size: 18ml

Intended use:

Storage: 2-8°C

Normal Mouse serum blocking buffer is used in immunohistochemical staining procedures to block non-specific staining when the host species of the secondary antibody is Mouse.

Kit components:

| Catalog No. | Size |
|---------------|-------------------------|
| NB-23-00187-1 | 100 mL (Ready to use) |
| NB-23-00187-2 | 18 mL (Ready to use) |

Recommended protocol:

1. Apply 100uL or enough volume of **Normal Mouse Serum Blocking Buffer** on tissue before adding primary antibody and incubate for 10 minutes at room temperature.
2. Blot off Normal Mouse Serum Blocking Buffer Solution (**Do not rinse**) and add primary antibody.

Precaution:

Please wear gloves and take other necessary precautions.

FOR RESEARCH USE ONLY

Neo-Biotech

74 rue des Suisses – 92000 Nanterre