

Neo Vero, Chemically Defined Medium, w/o L-Glutamine

#Cat: NB-58-0114 Size: 500ml

Product Information

Neo Vero

Chemically Defined Medium, w/o L-Glutamine, sterile-filtered

Cat. NO. NB-58-0114

General Information

Neo Vero is a chemically defined medium that promotes the growth of Vero cells for the production of viruses or recombinant proteins. Furthermore, Neo Vero is serum-free and animal/human origin-free.

Product Specifications

Appearance	Clear Red Solution
Specifications	Chemically Defined Serum-Free Animal Origin-Free Human Origin-Free
Storage and shelf life	Store at +2°C to +8°C protected from light. Once opened, store at +4° C and use within 6-8 weeks.
Shipping Conditions	Ambient
Concentration	1 X
Buffer system	NaHCO ₃ (2.2 g/L) for 5% CO ₂ Atmosphere
Formulated with	Epidermal Growth Factor (EGF), recombinant Insulin, recombinant Sodium Pyruvate Phenol Red Hypoxanthine Thymidine

For a greater stability and extended shelf life the medium is formulated without L-Glutamine. Antibiotics are not recommended; however, 10 ml/L of Antibiotic-Antimycotic (100X) containing penicillin, streptomycin, and amphotericin B may be used if required.

Instructions for Use

Adaptation:

For cells grown in serum supplemented medium or other serum-free medium little or no adaptation is needed and may be directly transferred to Neo Vero. It is advisable to keep a backup culture in the original media until cells have adapted. If suboptimal growth is observed after direct adaptation for 3–5 passages use the sequential adaptation method.

Sequential adaptation:

1. Subculture cells into a 25:75 ratio of supplemented Neo Vero to the original media. During the adaptation procedure seed at twice the normal seeding density ($2-8 \times 10^4$ viable cells/cm²).
2. Subculture cells when confluency reaches 70–90%. Subculture the cells in fresh pre-warmed 25:75 ratio

of supplemented Neo Vero to the original media. Once consistent cell growth with high viability has been achieved, passage cells into a 50:50 ratio of supplemented Neo Vero to original medium.

3. Repeat step 2 of this procedure, stepwise increasing the ratio of Neo Vero to original medium (75:25 followed by 90:10) until the cells are subcultured into 100% Neo Vero. Multiple passages at each step may be needed.
4. Continue to monitor and passage cells until consistent growth with high viability is achieved. After several passages in 100% Neo Vero, the culture is considered to be adapted.

Precautions and Disclaimer

This product is for research use only. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

Help Needed?

If you have any further questions regarding this product, please do not hesitate to contact our cell culture experts by email (info@neo-biotech.com) or phone (+33 9 77 40 09 09).